

EPview™ App - Release Notes

iOS

Version	Title	Description
3.7.8	Debugging	Stabilized camera connection. Faster App launch.
3.7.7	Debugging and updated About Text	Stabilized camera connection. Updated Copyright information and EULA.
3.7.2	Freehand Annotation	Implementation of Freehand Annotation. Further debugging.
3.6.8	Debugging and implementation of a calibration table	Stabilized the functionality of the QR code scanner. Implemented a calibration table to store up to 6 calibrations.
3.5.4	QR code scanner and updated the content of About Text	Stabilized the functionality of the QR code scanner. Updated About Text in the App: <ul style="list-style-type: none"> - Added a summary of the content - Added the notice of intended use
3.5.1	Launch	Launch of EPview App for iOS.

Android

Version	Title	Description
2.9.17	Debugging and usability improvements	Resolves an issue that made it difficult to add and remove control points to a polygon on some Android devices. Resolves an issue where the double tap gesture to enter a text annotation was not always immediately recognized on some Android devices. Text annotations are now automatically updated after the text has been entered. Fixed a bug that could cause the calibration dialogue to crash. Connecting to a camera no longer automatically activates the auto exposure mode. Fixed a bug that could cause issues with the exposure setting menu when activating the AC flicker reduction function.
2.9.10	Debugging and updated About Text	Updated Copyright information and EULA. Resolves an issue with the “redo” function not working properly. Resolves an issue where the App could crash if too high numbers were entered during length calibration. Updated QR code scanner. Added Android 13 support. Requires Android 8.1 or later.

2.9.6	Freehand Annotation	Implementation of Freehand Annotation. Stabilized the functionality of QR code scanner for Android.
2.8.4	Debugging and implementation of a calibration table	Stabilized the functionality of the QR code scanner. Implemented a calibration table to store up to 6 calibrations.
2.7.2	Updated the content of About Text	Updated About Text in the App: <ul style="list-style-type: none"> - Added a summary of the content - Added the notice of intended use
2.7.0	Launch	Launch of EPview App for android.

Windows OS

Version	Build	Title	Description
1.4	29191	Debugging and Windows 11 support	Resolves an issue where it was not possible in Ethernet LAN mode to manually connect to further EP50 cameras unless at least one camera was already found automatically by the software. Support of Windows 11
1.3	19857	Freehand Polyline	Implemented a freehand polyline to the drawings function.
1.2	19853	Debugging	Correction of calibration between live and snapshot.
1.1	19821	Launch	Launch of EPview for Windows operating system.

OSD

Version	Build	Title	Description
V4.0	20230523	Debugging	Resolves an issue which could cause acquired movies to stutter on playback. Resolves an issue where saved files after formatting the SD card were only shown when the camera was restarted.
V3.3	20190202	Launch	Launch of EP50.